


BAX® System Q7 Version 5.2.0.36505 Software Patch


Replaces v5.2.0.32948

To install the 5.2 patch


1. Ensure the current version of the BAX® System Q7 software installed on the computer is version 5.2.0.32948. Open the BAX® application. Go to Help> About. You should see the following:

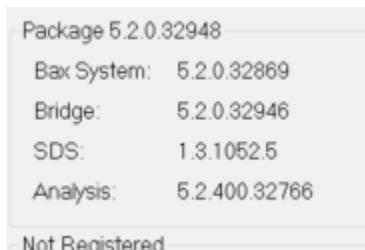

2. Close the application.
3. Reboot to ensure no Q7 components are active.
4. Download the zipped patch file (Q7_5.2_Update.zip) from the BAX System Q7 product page and copy to the following folder: C:\Program Files (x86)\BAX\Updates

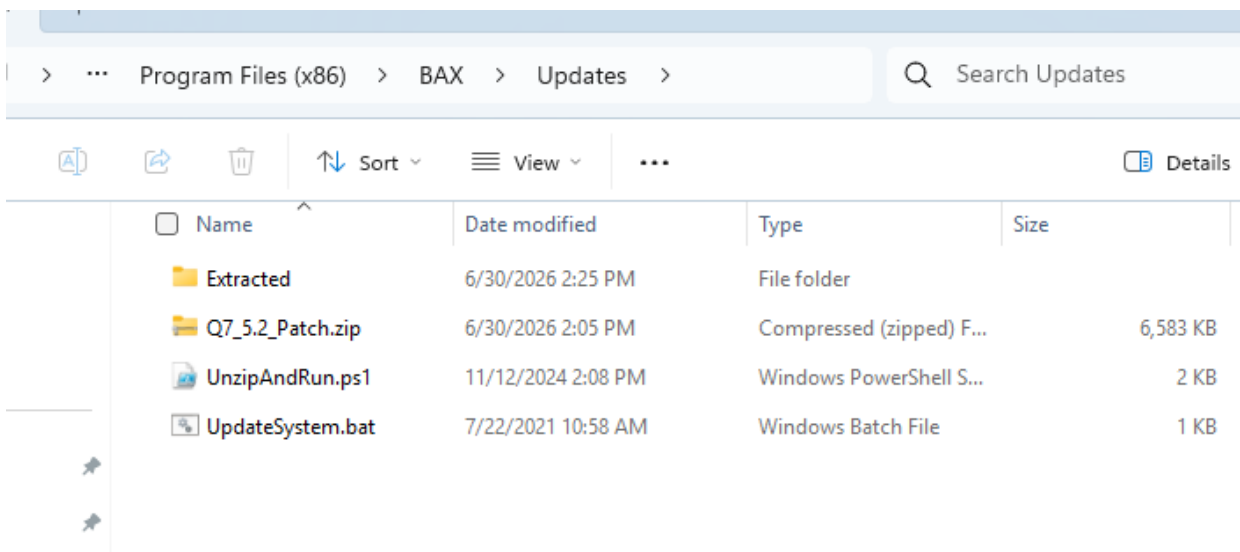

5. Execute UpdateSystem.bat by selecting: Run as administrator

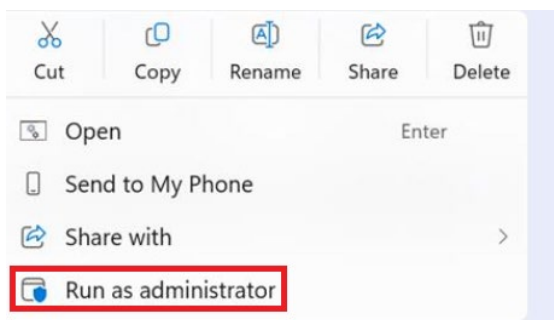

6. Execute the software: BaxSystem.exe –
 - a. See 5.2 hotfix version

Package 5.2.0.36505	
Bax System:	5.2.0.36504
Bridge:	5.2.0.36493
SDS:	1.3.1052.5
Analysis:	5.2.400.32766