



A rapid, qualitative, quantitative and reliable way to scan your GlutenTox® Sticks Plus for Reader

• Easy-to-use reader for standard lateral flow assays

• Provides quantitative, semi-quantitative & qualitative results

• Maintenance-free device

- Battery operated
- Small compact design
- Storage capacity of 100 scans
- Data transfer to a PC or laptop,
 with the Cube Data Reader software



How to use

1 Turn on device



To turn on the device press the button briefly, less than 1 second.

2 Ready for operation



The display shows ON and the reader is ready for operation.

The rapid test is now required.

3 Test insertion

Place the test into the adapter. If you use a strip adapter, you have to put the strip inside the adapter. If you use



a cassette adapter, you have to put the adapter on top of the cassette. Be sure that the adapter is correctly placed.

4 Measurement





Press the button briefly.

The display will show RFID or TEST.

5 Configuration data via RFID



Test configurations are uploaded via RFID before the measurement. Put the lot specific RFID card onto the top side of the device or hold the device against the RFID tag sticking on the test box until the configuration data has been uploaded (audible signal).



After the test is placed in the strip/cassette adapter, the reader can be placed on top. The slanted reader

corner and a corresponding counterpart inside the adapter help to find the correct position of the reader.



6 Measurement

Start the measurement by briefly pressing the button.



The device measures and the display shows RUN.



After a few seconds the result is displayed.



