

BAX® System Q7 Version 5.0.0.18610

Features and Instructions

BAX System Q7 software release v 5.0 contains several user interface features, new target options, and error logging improvements.

We strongly recommend that you install the latest version as soon as possible for optimal performance.

Please read all instructions thoroughly and ensure all required prerequisites are confirmed prior to beginning the installation of the software update.

If you currently have a Windows 7 computer system, please contact diagnostics.support@hygiena.com or your local Hygiena® representative for Windows 10 upgrade support.

This release is NOT applicable for Windows 7 or Windows 11 systems.

Software Assay Updates for Release v5.0.0.18610

- There are no changes to the analysis for BAX assays (vs. versions 4.21 and 4.22)
- Two foodproof® assays now with verified result interpretation (displayed as 'Certified' within the software):

| Product Name | Product Number |
|---|----------------|
| foodproof® <i>Aspergillus</i> Detection LyoKit (Real-time only) | KIT230145 |
| foodproof® <i>Salmonella</i> plus <i>Cronobacter</i> Detection LyoKit | KIT230131 |

foodproof assays with verified result interpretation do not require visual verification of the graphs. However, as a best practice, Hygiena recommends customers review all results.

- Seven additional foodproof assays with **experimental** analysis files for assistance with result interpretation:

| Product Name | Product Number |
|--|----------------|
| foodproof® <i>Listeria</i> Genus Detection LyoKit | KIT230088 |
| foodproof® <i>Listeria monocytogenes</i> Detection LyoKit | KIT230092 |
| foodproof® <i>Listeria</i> plus <i>L. monocytogenes</i> Detection LyoKit | KIT230129 |
| foodproof® <i>Salmonella</i> Detection LyoKit | KIT230099 |
| foodproof® <i>Salmonella</i> Enteritidis & Typhimurium Detection LyoKit | KIT230106 |
| foodproof® <i>Staphylococcus</i> Detection LyoKit | KIT230103 |
| foodproof® STEC Screening LyoKit | KIT230077 |


- foodproof/microproof® assays with improved **experimental** analysis files for assistance with result interpretation:

| Product Name | Product Number |
|--|----------------|
| foodproof® <i>Enterobacteriaceae</i> plus <i>Salmonella</i> Detection LyoKit | KIT230137 |
| microproof® <i>Legionella</i> Quantification LyoKit | KIT230119 |
| foodproof® <i>Salmonella</i> Genus plus Enteritidis & Typhimurium Detection LyoKit | KIT230134 |
| foodproof® <i>Vibrio</i> Detection LyoKit (Real-time only) | KIT230117 |
| foodproof® Yeast and Mold Quantification LyoKit | KIT230112 |

Important Note: All foodproof and microproof assays with **experimental** kit analysis files are designed to assist in result interpretation but require visual verification of the graphs.

Additional information for viewing data from **experimental** analysis files:


- Analysis files are now available for all assays (i.e., BAX, foodproof, microproof). By default, all assays with experimental analysis files will have analysis files set to “None”. The user can go to Operation → Analysis Settings to set the analysis file (Algorithm) to “Experimental” for targets that do not have verified analysis files. This setting enables improved assistance with results interpretation.

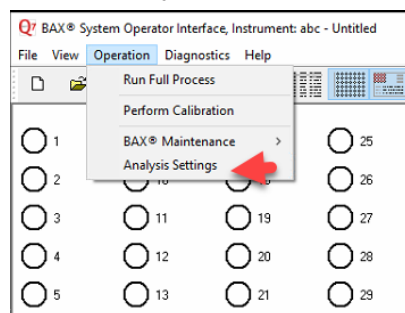


- All verified assays will have only one setting displayed as “Certified.”
- If you are using 4.22 to run foodproof assays, your selected setting will remain the same. Only the new foodproof assays in 5.0 will need to be adjusted to your preference.

Software Feature Updates for Release v5.0.0.18610

- Users are now able to register the Q7 to the SureTrend® application available on the cloud.
 - Ability to review data online both graphically and by result status with the SureTrend software.
 - To use this feature, you must register a SureTrend account <https://suretrend.hygiena.com/Register>
- Users can now export comma-delimited text files (.CSV)
- There is a one-step compressed log file export
- Software will now check that the correct calibration file is imported for the instrument in the case of equipment replacement by loaners or service units. Each Q7 has a unique calibration file and should only be used with that Q7.
 - Every calibration file generated with software version 4.0 and higher collects the serial number of the camera of the Q7 and the serial number of the thermal block. This information is stored in the calibration file. When a calibration file is imported, the application will compare the serial numbers in the calibration file to the serial numbers of the instrument on which you are importing. If the serial numbers don't match, the following pop-up message will appear. If this message appears, locate the correct calibration file (if it's a recent replacement unit, it can be found in the sleeve on the back of the unit - or - contact Hygiena to request the correct file). If the correct file cannot be located, calibrate the instrument as instructed in the BAX Q7 User Guide.

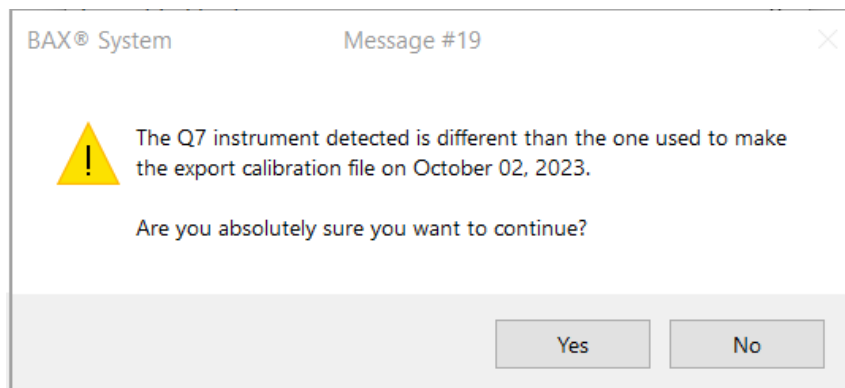

If "Yes" is clicked in the pop-up message above, a secondary message will appear

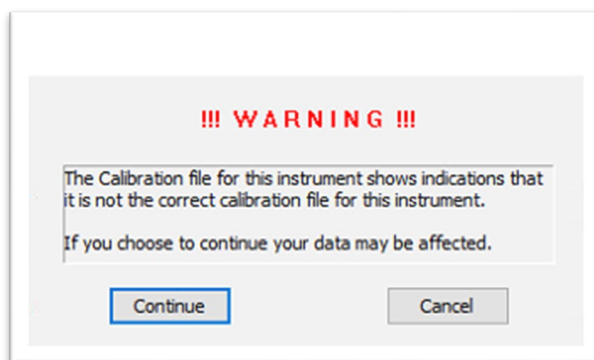

If 'Continue' is selected, the calibration will import, and no additional warnings or notifications will appear.


Software Upgrade Instructions:

1. Required Prerequisites:


- **Regional Settings are in English**
- **Core Isolation setting should be turned off in windows security settings**
- **Read/Write Permissions for the below file locations for all users**
 - C:\Program Files\MATLAB
 - C:\Program Files (x86)\BAX
 - C:\ProgramData\BAX Q7 System
 - All Power and Sleep Settings are set to Never.
- **Obtain HEX Calibration Kit (Product No.: KIT230340) from Hygiena if you plan to run foodproof/microproof assays that utilize the HEX dye (see kit instructions).**
 - This additional calibration plate is a stand-alone product and is not included with other calibration kits.
 - Please visit the [Contact Us](#) page for ordering and technical support.

2. Calibration storage and removal of current software:


- **Back up the system's calibration file.**
 - Select Operations > BAX Maintenance > Export Calibration from the menu bar.
 - Navigate to a new folder in which to store the backup file and click Save.
 - Name the calibration file (example: Q7Cal_date.bqc).
- **Uninstall the previous version of the BAX Q7 application.**
 - Navigate to Windows Settings and select Apps.

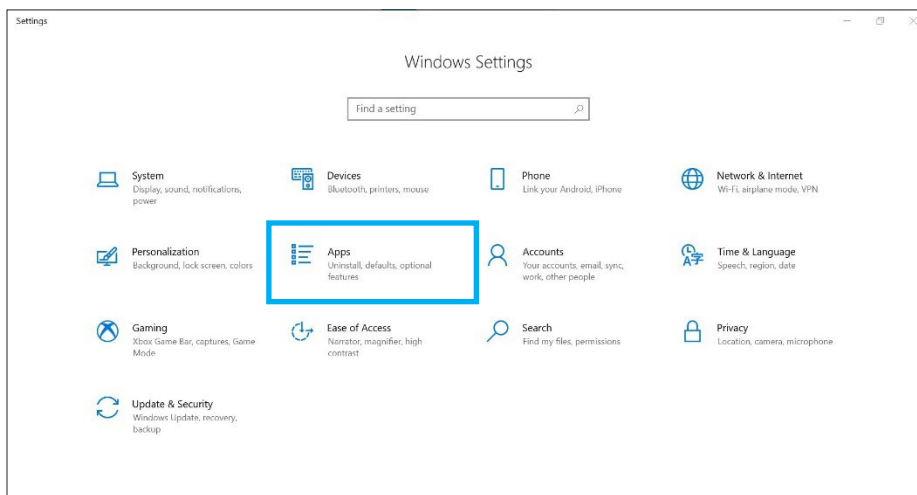

- Click on BAX System Q7, select Uninstall, and agree (click Yes) on the next window to initiate removal.

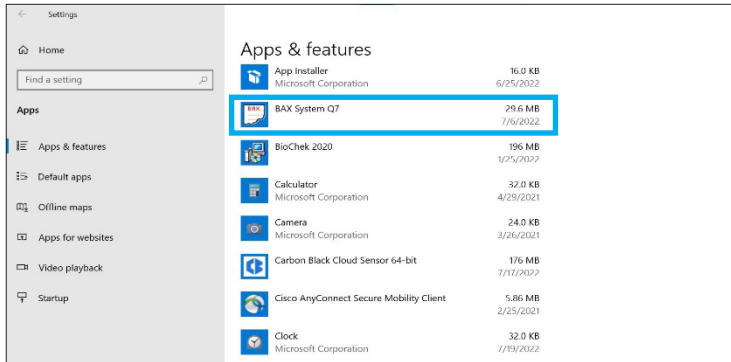


- If updating from 4.21 or 4.22, removal of MATLAB is not required. If updating from any previous versions, click on the MATLAB application, select Uninstall, and click Agree on the next window (this may take several minutes).

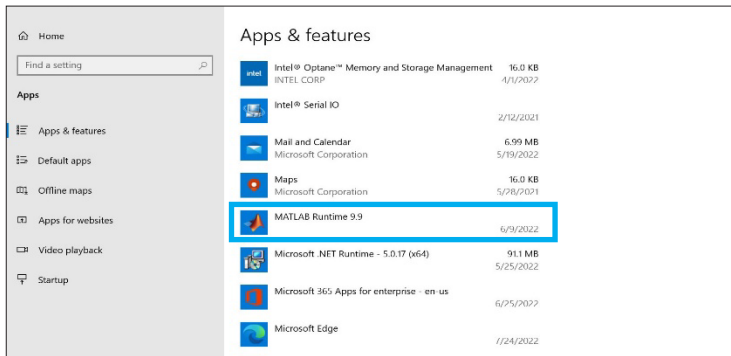


Restart computer **BEFORE** continuing to step 3.

3. Download 5.0.0.18610 software by clicking this [link](#) or the button from the BAX Q7 Page.

- Transfer download to a memory/USB stick if computer doesn't have internet access.
 - Click on the Q7 5.0.0.18610 icon and the program will automatically start execution of install.

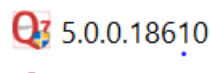


- A Windows protect screen may pop up. Select More Info, then select Run anyway.
 - On the next pop-up, select Yes to allow installation to begin

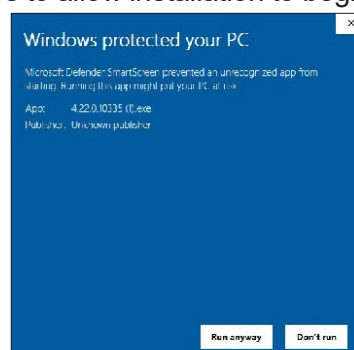

4. There may be some prerequisites that need to install to prepare your computer the BAX application installation.

If you see these, you may have to restart multiple times to get the prerequisites installed before MATLAB Runtime can install

BAX System Q7 - InstallShield Wizard

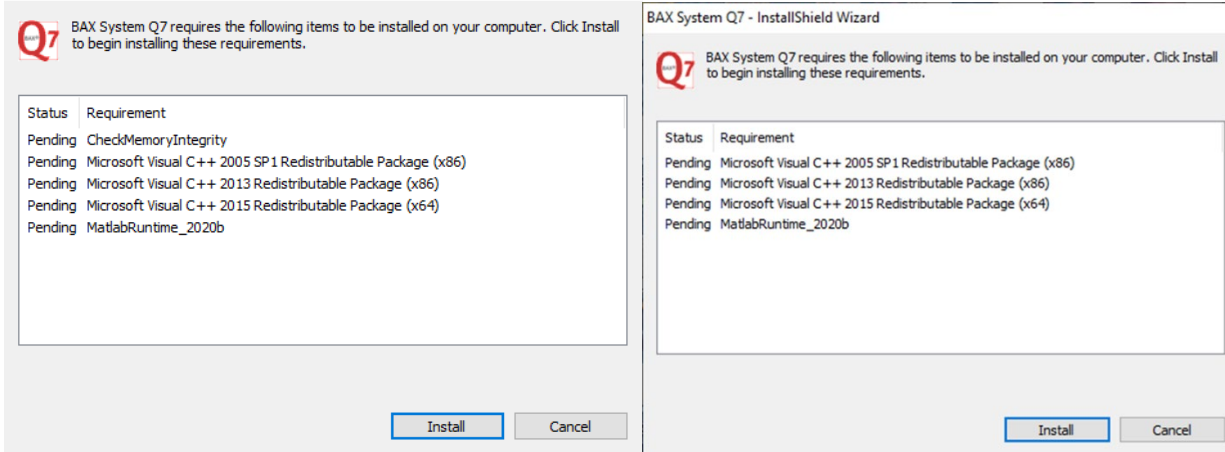

Checking memory integrity and Correct regional settings (not shown) are the first two prerequisites that need to be addressed or the Application will not install successfully.




In case this install error occurs: VCRUNTIME140.dll was not found (shown below).





- **Please download and install the patch: [vc_redist.x64.exe](https://www.microsoft.com/en-us/download/details.aspx?id=52685) from <https://www.microsoft.com/en-us/download/details.aspx?id=52685>. The installation will end when you click OK.**



5. **MATLAB updates will be installed (after prerequisites). Follow prompts below and select the default responses:**


After installing this patch, MATLAB Installation is continued.

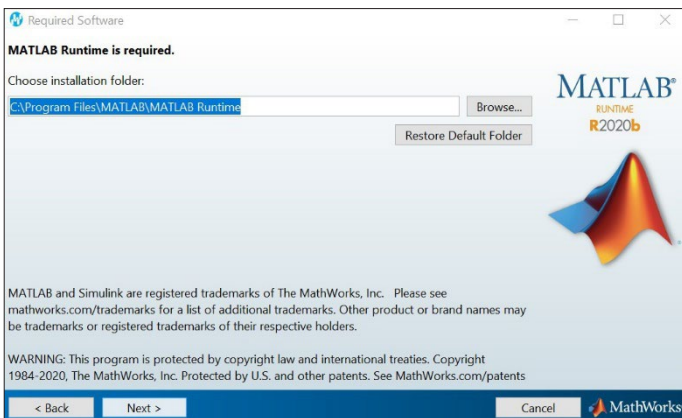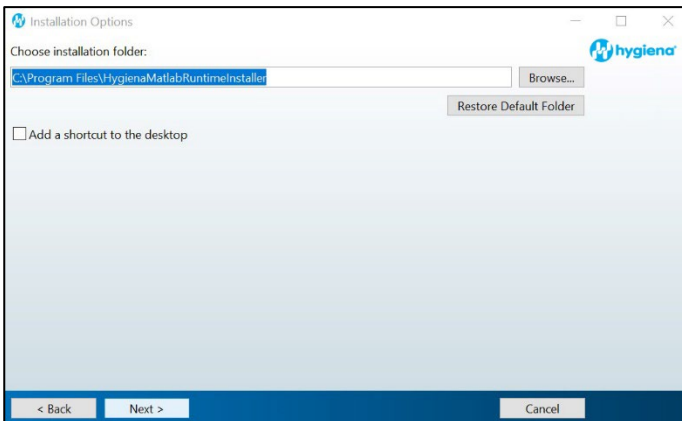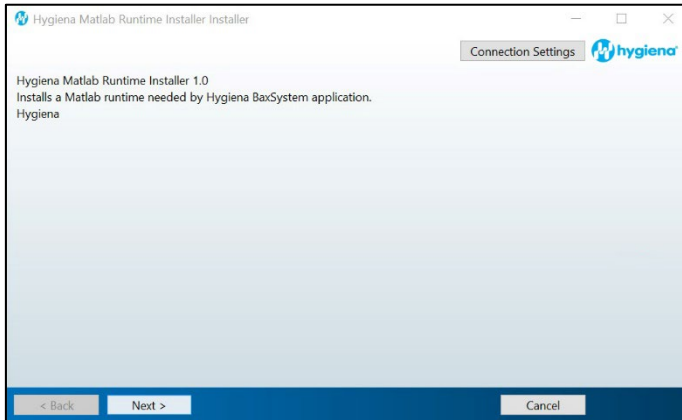

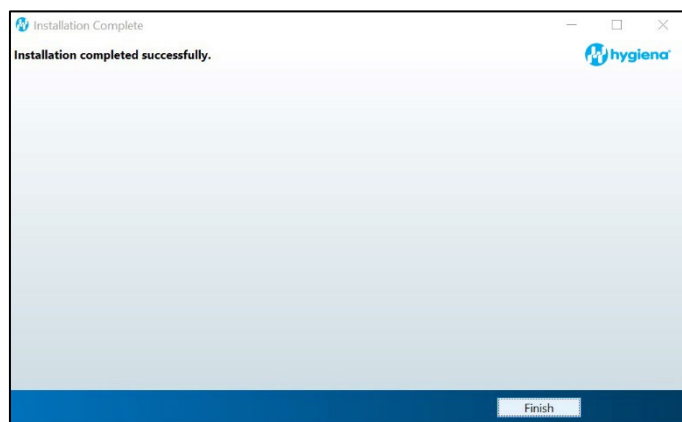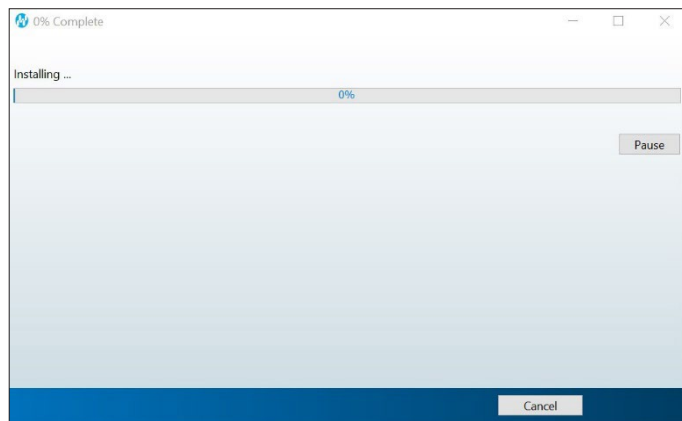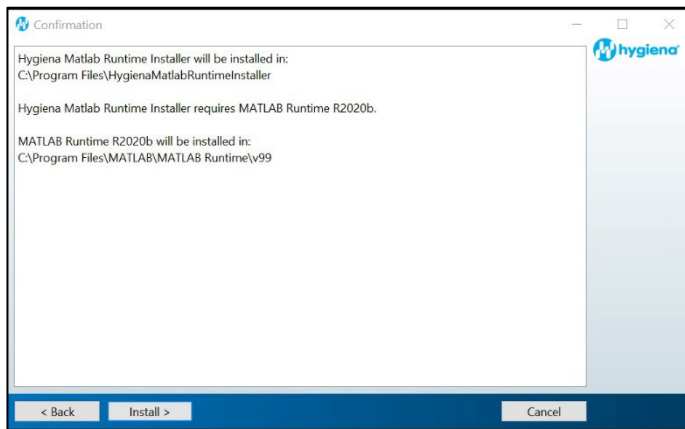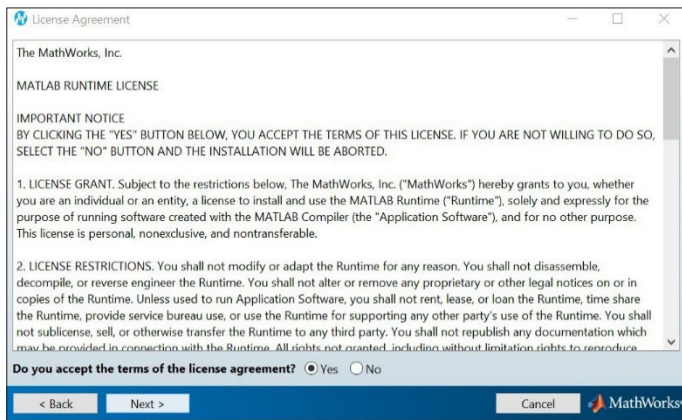


6. Following MATLAB installation, the BAX Q7 application installation will be initiated. Read the prompts carefully. Choose the default responses.

- Welcome Window – click Next.

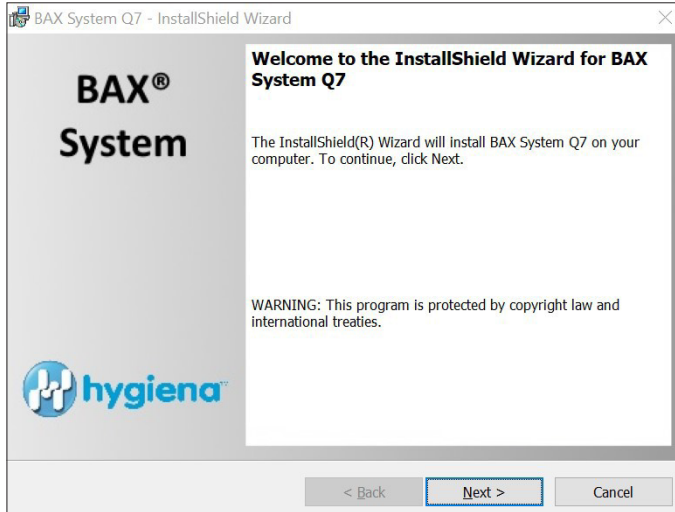


■ Calibration File Warning – ENSURE YOU HAVE BACKED UP THE CALIBRATION FILE.

- Check the box and click Next.

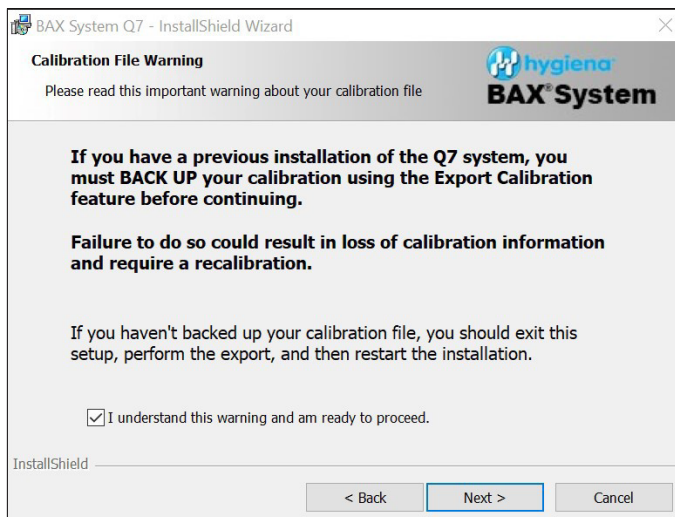


■ License Agreement.

- Check Accept Terms box and click Next.

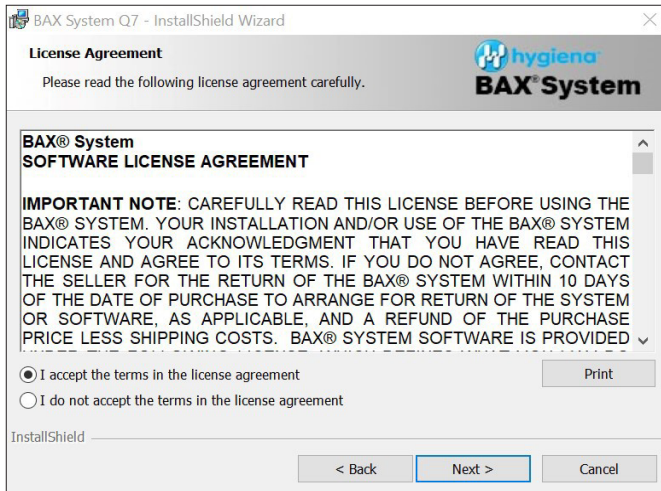

■ Custom Setup.

- Click Next.

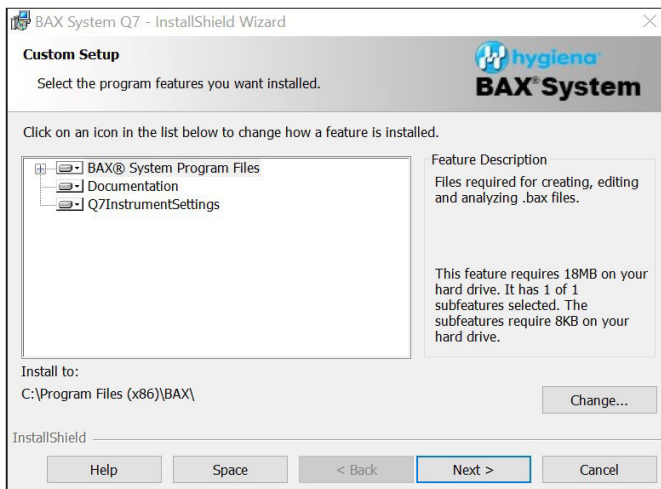


■ Password Protection.

- BAX default is disabled password protection. Choose an option for your lab, then click Next.





■ Ready to Install Program.

Click Install.

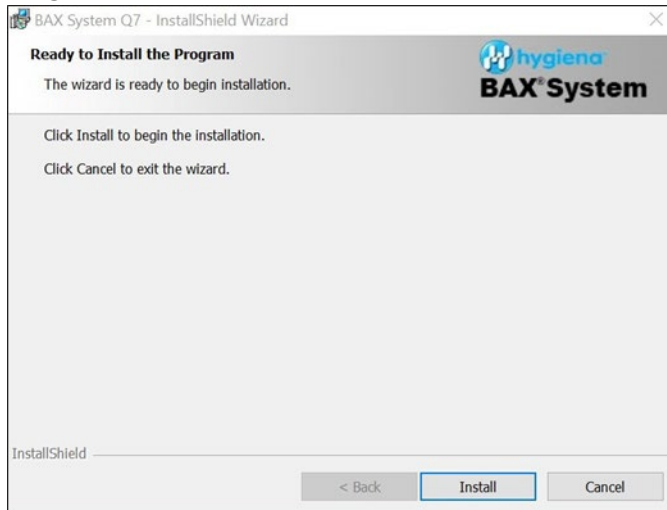


■ Installing BAX System Q7 Software.

- During installation, an Error Log Reporting window will pop up. We recommend NOT checking the box so the system can collect software and/or hardware errors in the Log Files to assist Diagnostics Support in troubleshooting. However, please review and decide what is appropriate for your lab.
- Select Next to complete the installation.

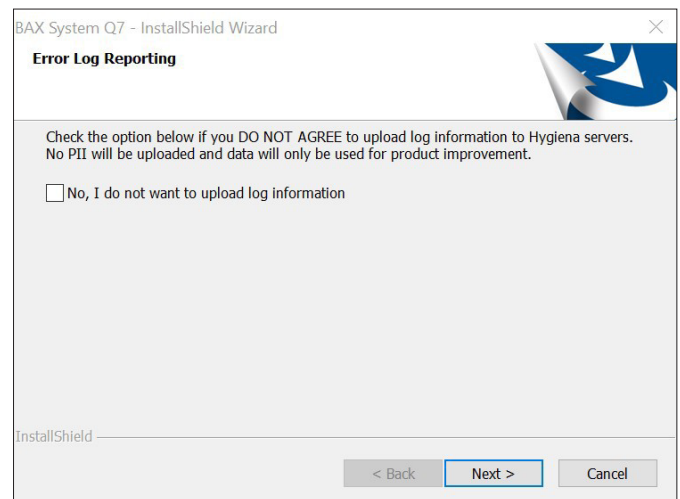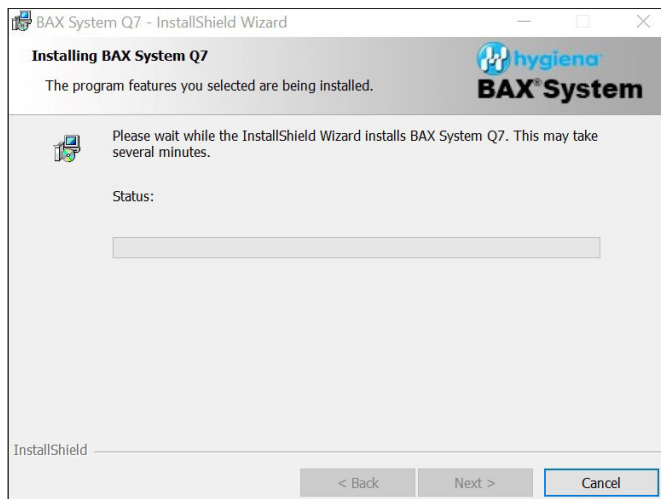


■ InstallShield Wizard Complete.


- Select Finish.

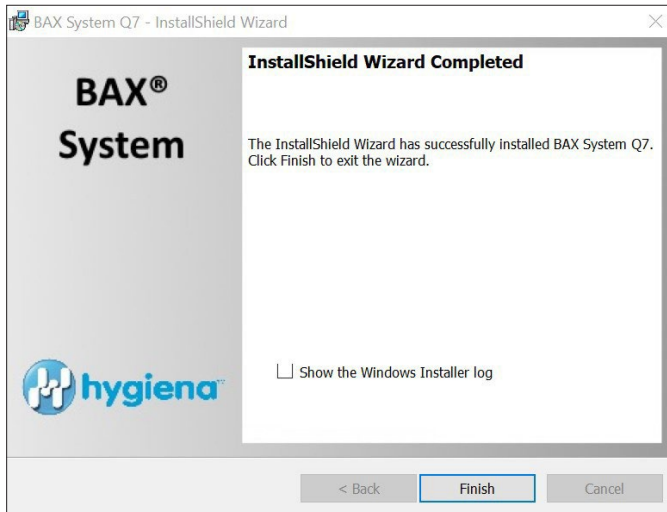

- Click on the desktop icon to initiate the program. The new open screen will have software version 5.0.0.18610.

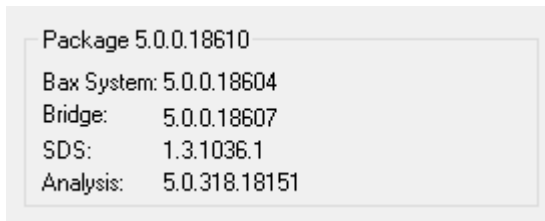

- Enter Instrument Name and Serial Number.
 - This will appear on reports and should be utilized.

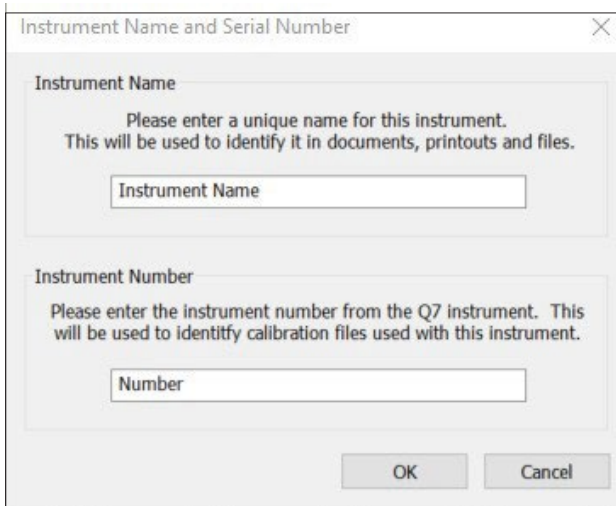

7. Check system's calibration


■ From the menu bar, select **Diagnostics > Calibration** to open the **Q7 Calibration Diagnostics** window

Click the Check Calibration button to determine the status:

- If calibration is available, no further actions are needed.
- If calibration is **NOT** optimal or **NOT** found:

Import the backed-up calibration file by selecting **Operations > BAX Maintenance > Import Calibration** from the menu bar and selecting the backed-up calibration file from step 2.

- To utilize many foodproof assays, additional calibration using the HEX calibration kit must be performed. Please follow the onboard instructions in the BAX Software to perform calibration.

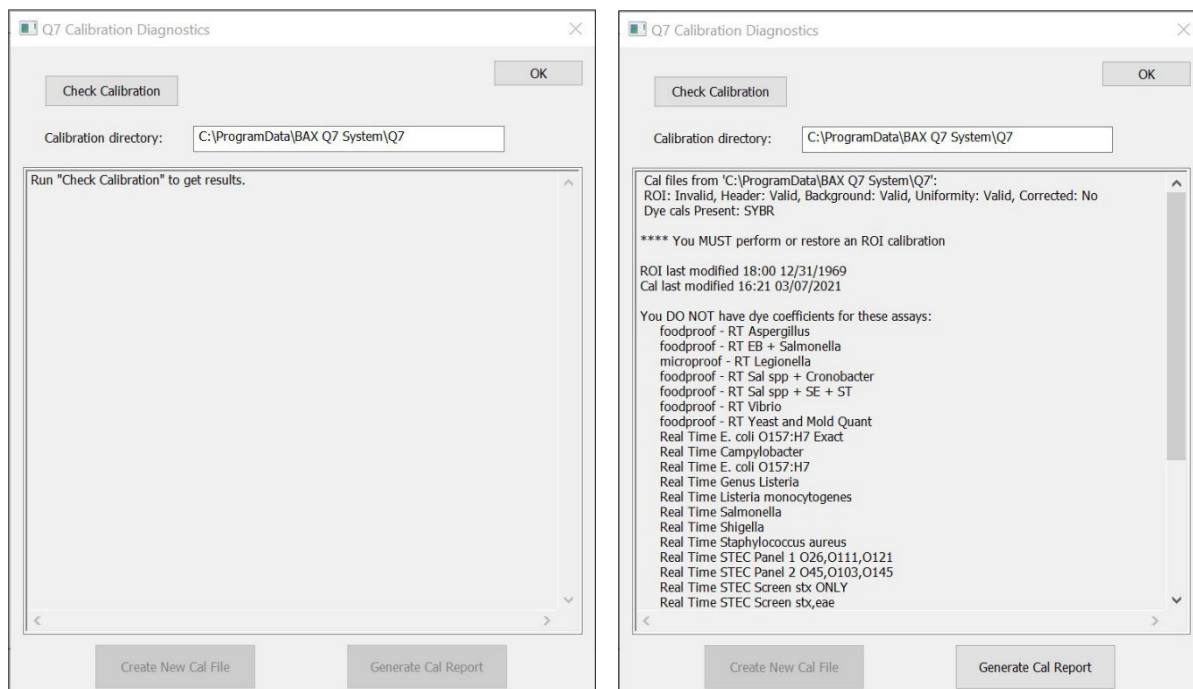

Appendix A: Algorithm Analysis across Software Versions

The table below provides information to identify algorithm consistency more easily across software versions and validations. There have been no changes to the assay-specific algorithm since the analysis version listed.

| Product Number | Assay | BAX Q7 software version with same algorithm as v5.0 |
|----------------|--|---|
| KIT2000 | BAX System Real-Time PCR Assay for <i>E. coli</i> O157:H7 | 4.0 |
| KIT2001 | BAX System PCR Assay for <i>Cronobacter</i> (<i>E. sakazakii</i>) | 4.0 |
| KIT2002 | BAX System PCR Assay for <i>L. monocytogenes</i> 24E | 4.0 |
| KIT2003 | BAX System PCR Assay for Genus <i>Listeria</i> 24E | 4.0 |
| KIT2004 | BAX System PCR Assay for <i>E. coli</i> O157:H7 MP | 4.0 |
| KIT2005 | BAX System Real-Time PCR Assay for <i>L. monocytogenes</i> | 4.0 |
| KIT2006 | BAX System Real-Time PCR Assay for <i>Salmonella</i> | 4.0 |
| KIT2007 | BAX System Real-Time PCR Assay for <i>Shigella</i> | 4.0 |
| KIT2008 | BAX System Real-Time PCR Assay for STEC Panel 1 | 4.0 |
| KIT2009 | BAX System Real-Time PCR Assay for STEC Panel 2 | 4.0 |
| KIT2010 | BAX System Real-Time PCR Assay for <i>Vibrio</i> | 4.0 |
| KIT2011 | BAX System PCR Assay for <i>Salmonella</i> 2 | 4.0 |
| KIT2016 | BAX System PCR Assay for Genus <i>Listeria</i> | 4.0 |
| KIT2017 | BAX System PCR Assay for <i>L. monocytogenes</i> | 4.0 |
| KIT2018 | BAX System Real-Time PCR Assay for <i>Campylobacter</i> | 4.0 |
| KIT2019 | BAX System Real-Time PCR Assay for Genus <i>Listeria</i> | 4.0 |
| KIT2020 | BAX System Real-Time PCR Assay for <i>Staphylococcus aureus</i> | 4.0 |
| KIT2021 | BAX System Real-Time PCR Assay for STEC Screening | 4.0 |
| KIT2039 | BAX System Real-Time PCR Assay for <i>E. coli</i> O157:H7 Exact | 4.0 |
| KIT230131 | foodproof <i>Salmonella</i> plus <i>Cronobacter</i> Detection LyoKit | 5.0 |
| KIT230145 | foodproof <i>Aspergillus</i> Detection LyoKit (Real-time only) | 5.0 |