



How to Install a BAX® System Q7 Software Patch

A patch is a rapid way to provide an easy-to-install solution for existing software located on the BAX® System Q7 computer. Patches must be overlaid on top of the proper existing software.

- Be sure to view the current software version within the Q7 software by navigating to the “Help” tab, then “About BAX System”.
- A pop-up window will appear, and the version can be viewed as the Package number.
- Utilize the package number to select the correct patch below.

4.21.3.9366 Patch

Instructions

1. Previous Q7 Install required (4.21.3.9366 version)
2. Make sure Q7 is not running by doing a reboot
3. Download the patch [HERE](#)
4. Copy patch (**4.21.3_CompletionCheck.zip**) to C:\Program Files (x86)\BAX\Updates
5. Execute C:\Program Files (x86)\BAX\Updates\UpdateSystem.bat as Administrator
6. Click on BAX System Q7 icon to run Q7

4.22.1.12282 Patch

Note: This hotfix includes the previous hotfix (4.22.1_CompletionCheck.zip)

Hotfix Patch Instructions:

1. Previous Q7 Install required (**4.22.1.12282 pkg or 4.22.010335 pkg**)
2. Make sure Q7 is not running by doing a reboot.
3. Download the patch [HERE](#)
4. Copy patch (5.0.1._MessagesUpdate.zip) to C:\Program Files (x86)\BAX\Updates
5. Execute C:\Program Files (x86)\BAX\Updates\UpdateSystem.bat as Admin
6. Click on BAX System Q7 icon to run Q7.

Versions Included:

Package	4.22.2.14335
Bax System:	4.22.1.12281
Bridge:	4.22.1.14333
SDS:	1.3.1036.1
Analysis:	4.22.0.9784

Differences between previous 4.22.1.12282 pkg:

1. 5944 - Create 2021.4 release patch/hotfix for the Q7 app not running.